



ROTOR BEATER MILL CR200

- Feed size □ ≤25mm
- Final fineness □ 40μm-10mm
- Speed □ 3000-10000rpm

The Rotor Beater Mill CR200 can grind both soft and hard materials through a combination of impact and shear actions, resulting in effective grinding.

APPLICATIONS

Pharmaceutical products, seeds, building materials, soil, chemicals, medicines, fertilizers, feed granules, grains, spices, coal, etc.

Rotor Beater Mill CR200

Working Principle

Samples enter the center of the grinding chamber through the hopper and are crushed by the rotor and sieve through a combination of impact and shear forces. When the sample size becomes smaller than the sieve's aperture size, it is directed into the receiving container. The mill door features a quick-locking design, providing easy access to the grinding chamber for convenient cleaning and rapid maintenance.

Rotor Beater Mill CR200

Accessories

Rotor Beater Mill CR200

180° bottom sieves

Suitable for soft, medium hard shell grain samples

Rotor Beater Mill CR200

360°bottom sieves

Suitable for soft and brittle grain samples

Rotor Beater Mill CR200

Rotor for CR200:

Mainly used for crushing grain, for example: wheat, corn, soybean.

FEATURES

- Adjustable speed ranging from 3,000 to 10,000 RPM to accommodate various application requirements.
- Maximum feed size of 25 mm.

- Achieves a final fineness of $d_{90} < 50 \mu\text{m}$.
- Equipped with a 180-degree sieve suitable for hard samples like grains with shells.
- Features a 360-degree sieve suitable for medium-hard, ductile, and fibrous samples.
- Provides a defined final fineness, thanks to sieves with aperture sizes ranging from 0.08 to 10.0 mm.
- A strong airflow generated by the high rotational speed rapidly cools the sample and aids in sample collection and cleaning of the grinding chamber.
- Easy cleaning facilitated by exchangeable push-fit grinding sieves and rotors.

Connect cyclone separator, small volume sample receiver

When Rotor Beater Mill CR200 is paired with a cyclone separator, it becomes well-suited for processing lightweight or small-volume samples. The inclusion of a cyclone separator broadens the Rotor Beater Mill's range of applications.

This enhancement not only efficiently addresses heat transfer issues during the grinding process, making cleaning a straightforward task, but also enhances the consistency and reproducibility of grinding outcomes.

Rotor Beater Mill CR200 Rotor Beater Mill CR200

TECHNICAL SPECIFICATIONS

Feed size	≤25mm
Final fineness	40μm-10mm
Speed	3000-10000rpm
Collector capacity	0.25-30L
Collector	0.25L□ 0.5L□ 1L□ 3L□ 5L□ 30L

Peripheral speed	21.4-71.2m/s
Sieves	0.08/0.20/0.25/0.50/1.00/2.00/4.0/6.0/8.0/10.0mm
Grinding chamber material	stainless steel, 1.1740 steel, hardened steel
Rated power	3.7KW
Power supply	380V□ 50/60Hz (110V Also available upon request)
Instrument size	590*705*1260mm
Package size	1000*1000*1450mm
Net weight	120kg



Torontech Inc.

251 Consumers Road, Suite 1200
Toronto, Ontario, M2J 4R3, Canada 🇨🇦

Phone: +1 416 368 2721 | Fax: +1 416 981 7652
Email: info@torontech.com

Torontech, LLC

601 Brickell Key Drive, Suite 700
Miami, FL 33131, USA 🇺🇸

Toll-Free: 1-866-383-7919
Email: sales@torontech.com

Torontech ME FZE

Meydan Grandstand, 6th Floor,
Meydan Road, Nad Al Sheba,
Dubai-261440, UAE 🇦🇪

Phone: +971 488 19252
Email: me@torontech.com

