



TORONTECH



AUTOMATED HOMOGENIZER - TORONAH™ 50

Automated Homogenizer - ToronAH™ 50

ToronAH™ 50 is an Automated Homogenizer designed to automate solvent addition, high-speed homogenization, and cutter/probe cleaning in one workflow. It replaces manual pre-treatment steps and supports unattended operation for high-throughput sample processing—up to 36 samples per batch.

ToronAH™ 50 is built for residue testing and routine lab prep where consistent homogenization and low carryover matter. Its flexible racks support multiple tube formats and compatibility with conventional centrifuge vials, helping standardize preparation across different methods and sample types.

APPLICATIONS

Automated Homogenizer - ToronAH™ 50 Applications

- Pesticide residue screening (food safety)
 - Automates homogenization and solvent addition for vegetable, fruit, and other agricultural samples used in multi-residue workflows.
- Veterinary drug residue testing
 - Supports consistent pre-treatment for animal-derived matrices such as meat, eggs, and milk, where repeatable homogenization is critical for extraction efficiency.
- High-moisture samples
 - Suitable for water-rich matrices like leeks, spinach, and celery, where uniform blending helps improve extraction repeatability.
- High-fat samples
 - Handles fatty matrices such as meat, sausages, and fish, supporting more stable preparation before downstream cleanup and analysis.


- High-fiber samples
 - Works well for fibrous materials including animal feeds and green forage, where strong homogenization is needed to reduce sample variability.




FEATURES

Automated Homogenizer - ToronAH™ 50 Key Features

- Fully automated workflow (low labor, high repeatability)
 - Automates solvent adding + homogenization + cutter cleaning, enabling unattended operation and reducing manual variability.
- High throughput for busy labs
 - Processes up to 36 samples per batch, improving daily sample capacity for routine screening.
- Cross-contamination control with automated probe cleaning
 - Generator probes can be cleaned automatically during operation, with multiple cleaning methods to help minimize carryover between samples.
- Flexible tube racks and broad compatibility
 - Multiple sample tube racks are available, and the unit is compatible with conventional centrifuge vials for easier integration into existing workflows.
- Safer, cleaner operation for solvent-based prep
 - Designed to save space in a fume hood and help reduce organic solvent exposure, with a transparent door and LED lighting for clear visibility.
- Built-in safety monitoring
 - Includes alarms for abnormal conditions such as cutter mis-position, motor overheat, and sample overload to support reliable daily operation.
- Easy graphical interface
 - Graphical software interface supports both experienced staff and new operators.

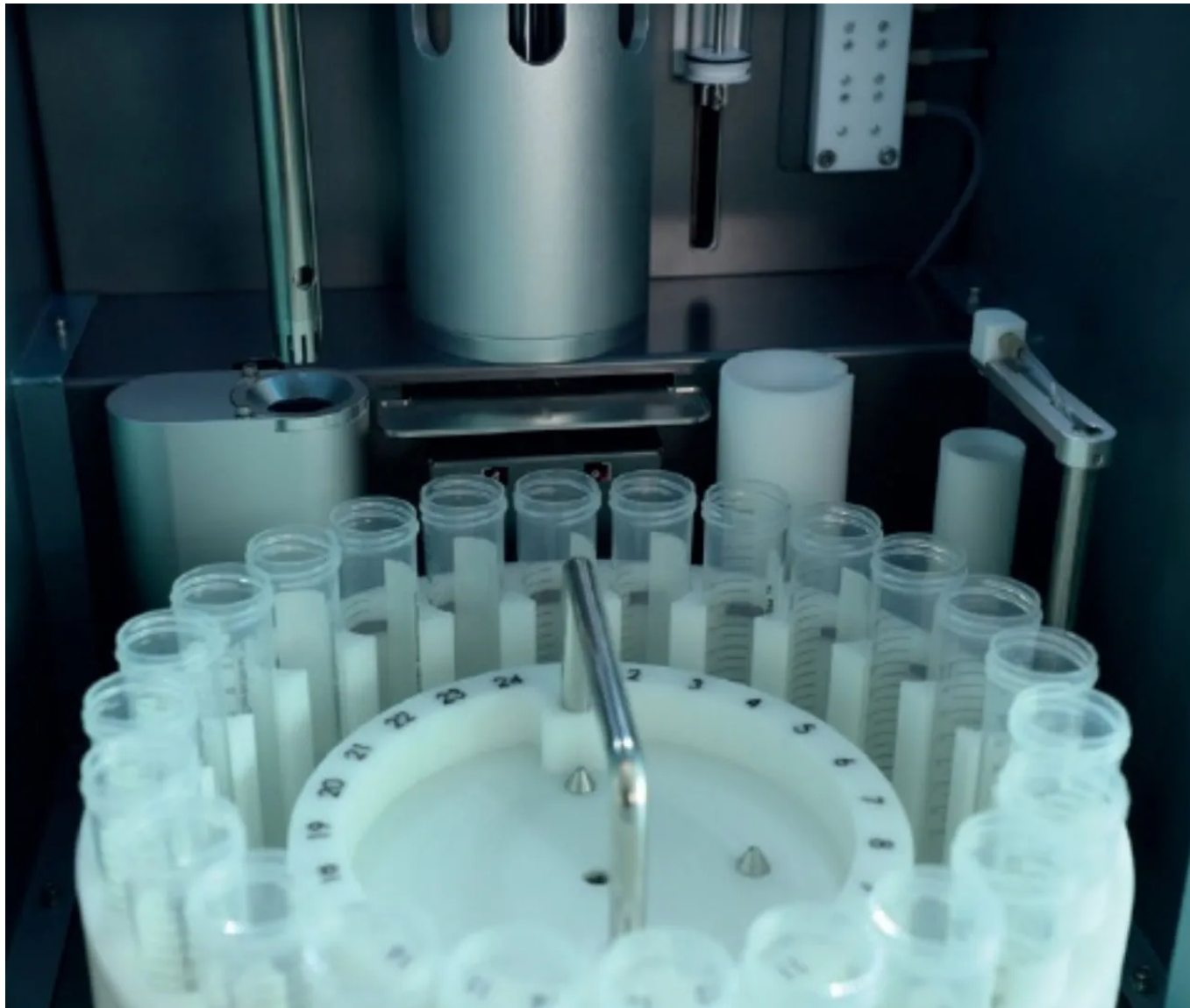

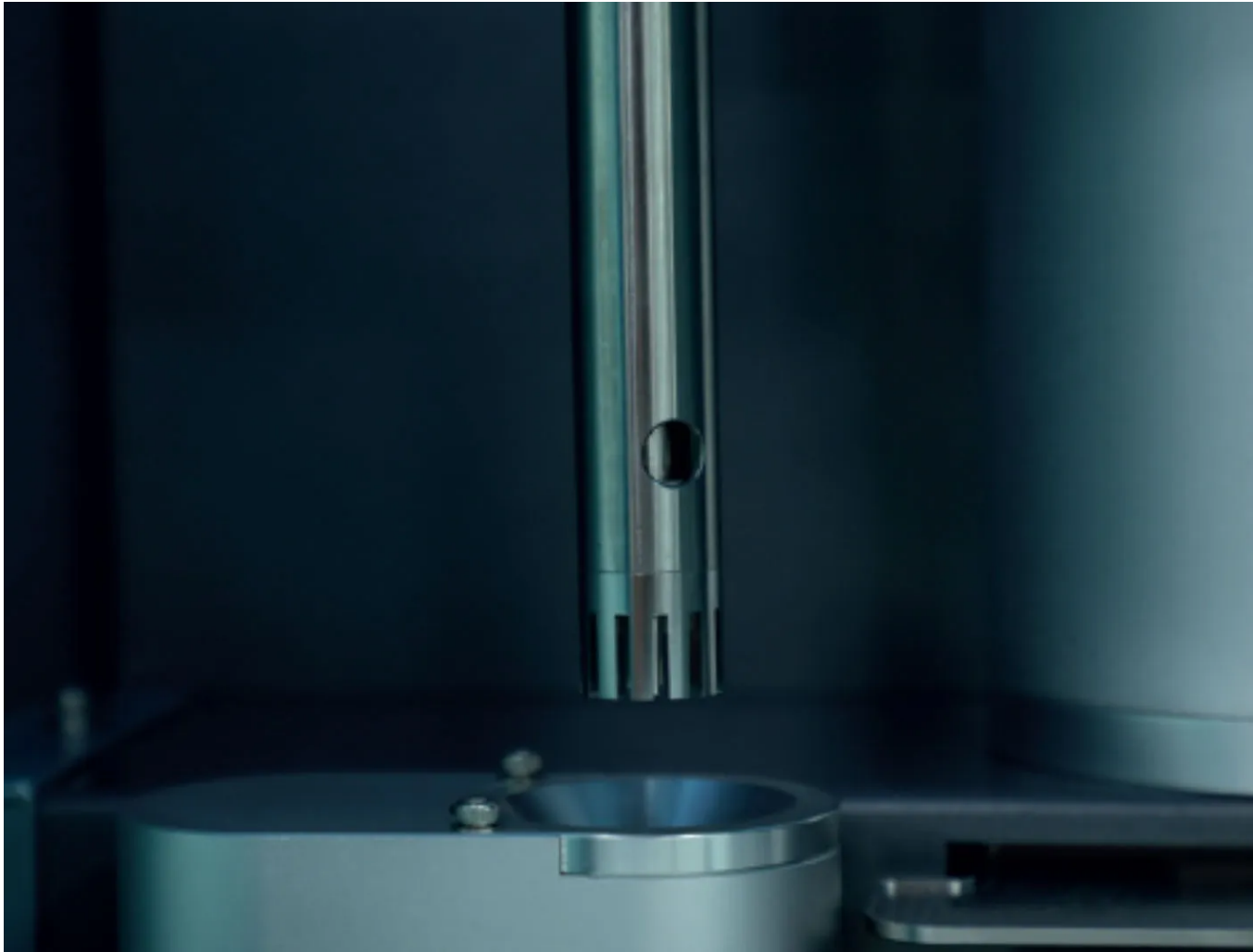

TECHNICAL SPECIFICATIONS

Automated Homogenizer - ToronAH™ 50 Technical Specification

Parameter	Specification
-----------	---------------

Carousel Rack Range	12, 16, 24, 36 positions; aluminum alloy, plastic
Oscillation Mode	Up-and-down vertical motion
Rinsing Mode	Water, organic solvents, ultrasound rinsing
Software Interface	10-inch built-in monitor screen control
Certificate	CE
Dimension	57 cm × 40 cm × 78 cm
Weight	75 kg
Power Rating	600 W
Speed Control	3,000 to 25,000 rpm
Generator Probes	10 mm × 108 mm; 19 mm × 204 mm
Sample Tube Range	15 mL, 50 mL, 100 mL, 120 mL, 200 mL



Torontech Inc.

251 Consumers Road, Suite 1200
Toronto, Ontario, M2J 4R3, Canada 

Phone: +1 416 368 2721 | Fax: +1 416 981 7652
Email: info@torontech.com

Torontech, LLC

601 Brickell Key Drive, Suite 700
Miami, FL 33131, USA 

Toll-Free: 1-866-383-7919
Email: sales@torontech.com

Torontech ME FZE

Meydan Grandstand, 6th Floor,
Meydan Road, Nad Al Sheba,
Dubai-261440, UAE 

Phone: +971 488 19252
Email: me@torontech.com

